BagMixer® 400

400 mL lab blenders



interscience

Why choose **BagMixer®** 400!

Best-seller

- Sold in more than 90 countries
- 35 years experience

High-quality

- Optimal analysis with the sterile BagFilter® and BagPage®
- In compliance with ISO 7218 standard
- Made in France by interscience

Brushless motor

- No wearing parts
- - 50% power consumption
- Extra silent

Safety

- The handle closes the blender and starts the blending
- Q-Tight[®]: no leaks from the top of the bag
- Security drip tray

BagMixer interscience

Easy cleaning

- All stainless steel
- Door opens at 270°
- Removable
 Click & Clean® paddles

PATENTED

Click & Clean® removable paddles

 Remove the paddles and easily clean the blending chamber

Guarantee

- Shock absorbers and window door life-time guarantee
- 3 year warranty (after registration)

PATENTED


Adjustable blending power

 Up to 28 kg of pressure: the hardest samples don't resist!
 e.g.: nuts, pills, seeds, dog biscuits etc.



Optimal extraction

- From as little as 30 seconds
- Easy control with the digital screen (speed, time, power)

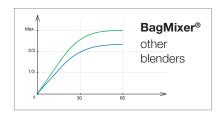

More than 15 000 BagMixer® in use every day!

Made in France by interscience,

the only manufacturer of blenders and sterile bags for microbiological analysis









WINDOW DOOR LIFETIME WARRANTY



BagMixer® 400 models:

The **BagMixer®** lab blender guarantees optimal extraction by the action of its two paddles. Adapted to any sample (food, agricultural, chemical, medical, environmental) to be analyzed between 50 mL and 400 mL, it guarantees homogenization in 30 to 60 seconds.









BagMixer® 400 S

Ref. 025 000

For more info on BagMixer® S & SW models, please refer to BagMixer® S & SW brochure.

> Silent blending



BagMixer® 400 P > The essential

Ref. 021 230

BagMixer® 400 W > Window door

Ref 022 230

BagMixer® 400 CC

> Click & Clean®

Ref 024 230

BagMixer® 400 SW

> Window door & LiquidSensor

Ref. 025 100

Also available: packs including bags and accessories; 100 mL and 3500 mL lab blenders. For more info, please contact us.

BagMixer® packs:



Starter Pack

- 1 BagMixer®
- + Sterile bags: 500 BagFilter® + 500 BagLight®
- + Accessories: 50 BagClip® + 1 BagOpen® + 1 BagRack®

Classic Pack

- 1 BagMixer®
- + Sterile bags: 500 BagFilter® + 500 BagLight® + 500 BagPage®
- + Accessories: 50 BagClip®
 - + 1 BagOpen® + 1 BagRack®
 - + 1 PetriPile® + 1 BagPipet®
 - + 1 000 BagTips® Regular
 - + 6 Clinet® 101

Premium Pack



- + Sterile bags: 1 000 BagFilter® + 2 000 BagLight® + 1 000 BagPage®
- ♣ Accessories: 50 BagClip®
 - + 1 BagOpen® + 2 BagRack®
 - + 2 PetriPile® + 1 BagPipet®
 - + 1 000 BagTips® Regular
 - + 6 Clinet® 101

BagSystem® complete range for sample preparation before analysis



BagFilter® Collect

Dilu**Flow**® Dilute



BagMixer® Blend



BagTools® Pipet

Technical specifications

		BagMixer® 400 P	BagMixer® 400 W	BagMixer® 400 CC
	Reference	021 230	022 230	024 230
	Removable and autoclavable paddles, Click & Clean® system	-	-	~
	Blending power indicator	-	-	✓
	Adjustable blending power (up to 28 kg of pressure)	-	-	~
S	Digital screen	-	-	✓
ture	Leak free Q-tight® system	-	-	✓
Key features	Side-by-side paddle stop	-	-	✓
Key	Silent brushless motor	-	-	✓
_	Window door	-	✓	✓
	Security drip tray	in option	✓	✓
	270° opening door	V	✓	✓
	All stainless steel	✓	✓	✓
	Compact and ergonomic	✓	✓	✓
SL	Door weight	1050 g 950 g		50 g
atio	Blending time	30, 60, 90, 120, 150, 180, 210 sec., ∞ from 1 sec. to 59 m		from 1 sec. to 59 min, ∞
Technical specifications	Blending speed	fixed speed: 8 strokes/sec.		variable speed: 4, 6, 8, 10 strokes/sec.
l sp	Dimensions (w x d x h)	27 x 38.5 x 28 cm		27 x 41.5 x 26.5 cm
ica	Weight	16.6 kg		16 kg
chr	Power	110-127/220-240 V~ 50-60Hz		100-240 V~ 50-60Hz
Pe	Capacity	50-400 mL		
es	Made in France	✓	✓	✓
Warranties	Life-time warranty	Unbreakable shock absorbers and window door		
	3 year warrabty (after registration)	✓	~	✓
Manufactured in compliance with		9001 2002/96/EC 2	RoHS 2002/95/EC MADE IN IRRANCE	

Discover our complete range for microbiology



BagFilter® Sterile bags



Dilu**Flow**®
Gravimetric dilutors



FlexiPump® Peristaltic pumps



easy**Spiral** Dilute® Automatic diluter et platers



Automatic colony counters & inhibition zone readers

Your local distributor

interscience

30, chemin du Bois des Arpents - 78860 St Nom - FRANCE T: +33 (0)1 34 62 62 61 - F: +33 (0)1 34 62 43 03 - Email: info@interscience.com

interscience usa & canada

3 2 Cummings Park - Woburn, MA 0 1 8 0 1 - USA T: +1 781 937 0007 - F: +1 781 937 0017 - Email: sales.usa@interscience.com

interscience CHINA

上海市徐汇区吴兴路277号锦都大厦588室 - 200030 电话: +86 (0)21-64739390 - +86 189 3097 0733 - 邮址: sales.china@interscience.com

interscience south-EAST ASIA

541 Orchard road - 09-01 Liat Towers - SINGAPORE 238881 T: +65 6933 1389 - +65 8549 1217 - Email: sales.asia@interscience.com



BagMixer® S&SW

400 mL lab blenders



Super Silent Blending!

interscience



Our quality for your lab

- Full range of lab instruments for microbiological analyses
- Designed and made in France
- Leadership in R&D for innovative products
- Worldwide distribution network in more than 85 countries

20 years of innovations for great sample preparation

With over **15,000 BagMixer**® sold around the world, we are recognized today as the world specialist for sample preparation with the **BagMixer**® lab blenders range and the **BagSystem**® range of products.

		1995	BagMixer® released World's 1st stainless steel window door lab blender	
		1998	MiniMix® 100 mL lab blender released Adjustable Blending Power to adjust the blending chamber to the sample 3 year warranty worldwide	TENTED
_ •	ar codes	1999	JumboMix® 3500 mL lab blender released Lifetime warranty on shock absorbers Lifetime warranty on window door	
_ •	Indiana.	2004	BagMixer® VW 400 mL lab blender released User-friendly features with the digital multifunction screen	
_ •	Beyldinor - 1 - ander 70 - 1 - ander 70	2007	Click & Clean® paddles Easy cleaning of the blending chamber with the removable paddles PA	TENTED
		2010	BagMixer® is a best-seller in more than 80 countries.	
-	Barg VI xox	2013	BagMixer® S and BagMixer® SW 400 mL lab blender released LiquidSensor alarm & TotalAccess System patents PA	ATENTED

Unique technology for an optimal bacterial extraction

Adjustable blending power -

Up to 28 kg of pressure per paddle

PATENTED



Small/hard samples: Move the paddles forward for more pressure



Big/soft samples: Move the paddles backwards for less pressure

Stainless steel body

High quality 304 L stainless steel

Robust and reliable

3 year warranty

Lifetime warranty on window

Lifetime warranty on shock absorbers

Advanced lighting system

LightCode System: indicates the state of the blending process (under the unit)

LED lighting: lights up the blending chamber during blending process and cleaning*

BagMixer® S



interscience

Ultra stable

Paulstra[™] anti-vibration feet Non marking

Super Silent Blending

Less than 48 dB

Noise cancelling casing

Double layer window door











SHOCK ABSORBERS LIFETIME WARRANTY WINDOW LIFETIME WARRANTY

Advanced new features



TotalAccess System


The cleaning of a blending chamber used to be time consuming. With the TotalAccess System, cleaning is now easy with the opening of the door to a completely flat position. Remove the Click & Clean® paddles and the drip tray: the bright polycarbonate chamber is easy to wipe thanks to its round corners.

PATENTED

Super Silent

Classic paddle blenders are very loud and induce fatigue. The BagMixer®S & SW will really change the work comfort of your lab. They are so silent that you can make a phone call next to them while blending: less than 48 dB in comparison with 54 dB with their closest competitor.

Max speed with 250 mL water + 22 g cheese. Tested by Laboratoire National d'Essai. Data available on request.

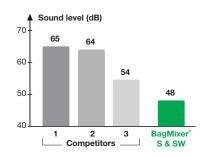



ISO 3744 acoustic test



LightCode System

Our new BagMixer® is so quiet that you won't even notice when the blending is over! We added some LED lights that illuminate the lab bench so that you know when the blending process is finished, even at a distance.

The light indicates the status:



Green: Ready







Red: LiquidSensor has detected a leak

LiquidSensor

If a blender bag leaks, the BagMixer® stops automatically. The sensor detects any liquid and alerts you right away, allowing you to save your sample and enjoy a quicker clean up!

PATENTED





EnergySaver

The new brushless motor technology allows the BagMixer® to consume 50% less energy than the brushed motor competitors. There are no wearing parts, so it can last much longer. The new AutoSleep function puts the BagMixer® in sleep mode after 30 minutes of inactivity.

BagMixer® 400 range

For the analyses of solid samples



BagMixer® 400 P

> The essential

Ref. 021 230



BagMixer® 400 W

> Window door

Ref. 022 230



BagMixer® 400 CC

> Click & Clean®

Ref. 024 230



BagMixer® 400 S

> Silent blending

Ref. 025 000



BagMixer® 400 SW

> Window door & LiquidSensor

Ref. 025 100

Also available: packs including bags and accessories; 100 mL and 3500 mL lab blenders. For more info, please contact us.

Complementary products & accessories





BagFilter® 400
Sterile bag with lateral filter

Ref. 111 425



BagPage® 400 Sterile bag with full-page filter Ref. 122 025



BagLight® 400 Sterile bag Ref. 132 025



BagSeal® Thermal sealing unit

Ref. 261 000



BagOpen® 400 Bag opener/holder

Ref. 211 040



BagRack® 400 & DripTray® (in option) Storage rack for bags Ref. 221 040



BagClip® 400
Closing clips

Ref. 231 040



BagPipet® & BagTips®

Pipet - **Ref. 251 091** Sterile tips - **Ref. 252 019**

3576 Bagkwars&S.W.B. 0216. Pictures and informations in this brochure are not contractually binding. Diul-Bow, BagFliter, BagLight*, BagFliter, BagTools* are trademarks. This document is the translation of the french brochure and is not contractually binding. INTERSCIENCE reserves the right to change or improve the specifications of its products without notice. Please visit www.interscience.if for current updates and additional information. RCS 950 356 220 Versailles. INTERSCIENCE SARL F78860.

Technical specifications

	BagMixer® 400 S	BagMixer® 400 SW	
Reference	025 000	025 100	
Adjustable blending power	✓	✓	
Side-by-side paddle stop	✓	✓	
Click and Clean® System: removable paddles	✓	✓	
304 L stainless steel body	✓	✓	
Integrated security stainless steel drip tray	✓	✓	
180 Watts brushless motor	✓	✓	
Multifunction digital screen	✓	✓	
4 adjustable speeds	stable speeds		
TotalAccess System: door opens flat	✓	✓	
LightCode: color status LED	✓	✓	
White polycarbonate blending chamber	✓	✓	
White LED light in blending chamber	olending chamber ✓		
Double layer window door	-	✓	
LiquidSensor: detects leaking bags	-	✓	
Sound level with 250 mL water + 22 g cheese	48 dB		
Door weight	1050 g	950 g	
Blending time	from 1 s to 59 mn, ∞		
Blending speed	4, 6, 8, 10 strokes/s		
Dimensions (w x d x h)	27.5 x 46.5 x 36 cm		
Weight	25.2 kg	25 kg	
Power	100-240 V∼ 50-60 Hz		
Capacity	from 50 to 400 mL		
Made in France	✓	✓	
3 year warranty (after registration)	✓	✓	
Lifetime warranty on shock absorbers	✓	✓	
Life-time warranty on window	-	✓	

 $\textbf{BagMixer}^{\texttt{o}} \text{ is delivered with: Power cord, user's manual, free sterile packs of filter bags.}$

Certified











Your local distributor

interscience INTERNATIONAL

30, chemin du Bois des Arpents - 78860 St Nom - FRANCE T: +33 (0)1 34 62 62 61 - F: +33 (0)1 34 62 43 03 - Email: info@interscience.com

interscience USA & CANADA

 $3.2\,$ C u m m i n g s P a r k - W o b u r n , M A 0 1 8 0 1 - U S A T: +1 781 937 0007 - F: +1 781 937 017 - Email: sales.usa@interscience.com

interscience CHINA

上海市徐汇区吴兴路277号锦都大厦588室 - 200030 电话: +86 (0)21-64739390 - +86 189 3097 0733 - 邮址: sales.china@interscience.com

interscience SOUTH-EAST ASIA

541 Orchard road - 09-01 Liat Towers - SINGAPORE 238881 T: +65 6933 1389 - +65 8549 1217 - Email: sales.asia@interscience.com





JumboMix® 3500

Large capacity lab blender



Accuracy and heavy duty blending for large samples

interscience

Blend up to 3500 mL!

JumboMix® lab blender guarantees optimal bacterial extraction by the action of its 2 paddles. Adapted to any volume between 200 mL and 3500 mL, it guarantees the homogenization in 30 to 60 s of all samples (up to 400 g).

Applications: *E. Coli O157:H7*, trichinella (WarmMix® model), *Salmonella spp*, baby milk powder, chocolate, sterility control tests, Cryptosporidium and Giardia parasites, sample pooling (e.g.: 15 samples of 25 g to reach 375 g)...

Powerful and heavy duty

- Up to 3500 mL capacity
- Ideal for your 375 g samples diluted to 1/10th
- Very silent

Ergonomic

- Side-by-side paddle stop: easy insertion of the bag
- All 304L stainless steel trolley, ergonomic height (in option)
- Digital screen

PATENTED

Easy cleaning

- All stainless steel
- 270° fully opening door
- Click & Clean® removable paddles

Security

- The handle shuts and starts the blending
- Q-Tight® closing: no leaks upwards
- Security drip tray



Guarantee

- Life-time warranty: Shock absorbers and window
- 3 year warranty (after validation of the card)
- Made in France

PATENTED

Adjustable blending power

- Adapt the blending power to your sample
- Up to 22.6 kg thrust: hard samples are blended




Before

After

Ideal for your applications

- 375 g for Salmonella spp
- 75, 325 or 375 g sample for *E. Coli O157:H7*
- 125 g sample for Listeria and Listeria monocytogenes
- Sample pooling

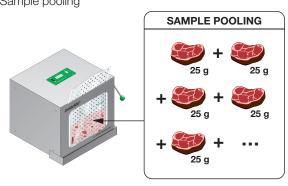

JumboMix® 3500

3 models available:



JumboMix® 3500 P CC

> Essential features

Ref. 031 230



JumboMix® 3500 W CC

> Window door

Ref. 032 230



JumboMix® 3500 WarmMix® CC

> Thermal regulation up to + 50°C

Ref. 033 230

ISO 7218









Technical specifications at the back of the brochure -

JumboLine®

Ideal for your preparation of 375 g samples diluted to 1/10th!





BagFilter®

Lateral filter bags for lab blenders

2000 mL - Ref. 111 200

3500 mL - Ref. 113 510

BagPage®

Full-page filter bags for lab blenders

2000 mL - Ref. 122 200

3500 mL - Ref. 123 010

BagLight®

Bags for lab blenders

2000 mL - Ref. 132 200

3500 mL - Ref. 133 025



JumboStand Ergonomic trolley

Ref. 037 000



BagOpen® 3500 Bag opener

Ref. 211 350



BagRack® 3500 Storage rack for bags

2 comp. - **Ref. 221 352** 4 comp. - **Ref. 221 350**



BagClip® 3500 Closing clips

Ref. 231 350



BagPipet® & BagTips®

Pipet - **Ref. 251 091**

Tips - Ref. 252 024

Technical specifications

		JumboMix® 3500 P CC	JumboMix® 3500 W CC	JumboMix® 3500 WarmMix® CC	
	Reference	031 230	032 230	033 230	
	Window door	-	✓	-	
	270° opening door	✓	✓	✓	
	Security drip tray	in option	✓	✓	
	Adjustable blending power (up to 22.6 kg thrust)	✓	✓	✓	
F	Blending power indicator	✓	✓	✓	
EQUIPMENT	Digital screen	✓	✓	✓	
	Q-Tight® anti-leak system	✓	✓	✓	
	Side-by-side paddle stop	✓	✓	✓	
	Click & Clean® removable paddles (patented)	✓	✓	✓	
	All 304 L stainless steel	✓	✓	✓	
	Compact and ergonomic	✓	✓	✓	
	Regulated heating door	-	-	✓	
	Door weight	6.9 kg	3.9 kg	6.1 kg	
¬ Si	Variable blending time	from 1 s to 59 min, ∞, count-down			
ĕ ĕ	Variable speed	1.5 / 3 / 4.5 / 6 strokes/s			
TECHNICAL SPECIFICATIONS	Dimensions (w x d x h)	52 x 47 x 47 cm		63 x 50 x 47 cm	
교등	Weight	53.2 kg	50.2 kg	63 kg	
SPE	Power	100-127 V~ / 220-240 V~ 50-60 Hz			
	Capacity	200 - 3500 mL			
Ţ	Made in France	✓	✓	✓	
3AN	Life guarantee	Shock absorbers and window door			
	3 years warranty (after registration)	✓	✓	✓	
Certified production		Products made for INTERSC by Interlab, an ISO 9001 certif	**	RoHS 2002/95/EC MADE IN FRANCE	

Delivered with: Power cable, user's manual, security DripTray*, packs of bags. (*only on JumboMix® 3500 W CC and JumboMix® 3500 W armMix® CC).

Discover our complete range for microbiology



insta**BAG**®
Blender bags
with dehydrated media



Dilu**Flow®**Gravimetric dilutors



FlexiPump® Peristaltic pumps



easy**Spiral** Dilute® Automatic diluter & platers



ScanStation®
Real-time incubator
& colony counter

Your local distributor

interscience

30, chemin du Bois des Arpents - 78860 St Nom - FRANCE T: +33 (0)1 34 62 62 61 - Email: info@interscience.com

interscience USA & CANADA

3 2 Cummings Park - Woburn, MA 0 1 8 0 1 - USA P: +1 781 937 0007 - F: +1 781 937 0017 - Email: sales.usa@interscience.com

interscience CHINE

上海市徐汇区肇嘉浜路798号坤阳大厦1903室- 200030 电话: +86 (0)21-64739390 - +86 189 3097 0733 - 邮址: sales.china@interscience.com

interscience ASIE DU SUD-EST

The Centropod, 80 Changi Rd - 05-07 Singapore 419715 T: +65 6909 0825 - M: +65 8118 5178 - E-mail: sales.asia@interscience.com