




Microbiology Focus

Volume 1.2, 2009



New Fast and Specific Detection Method of Pathogenic Fungi in Tissue

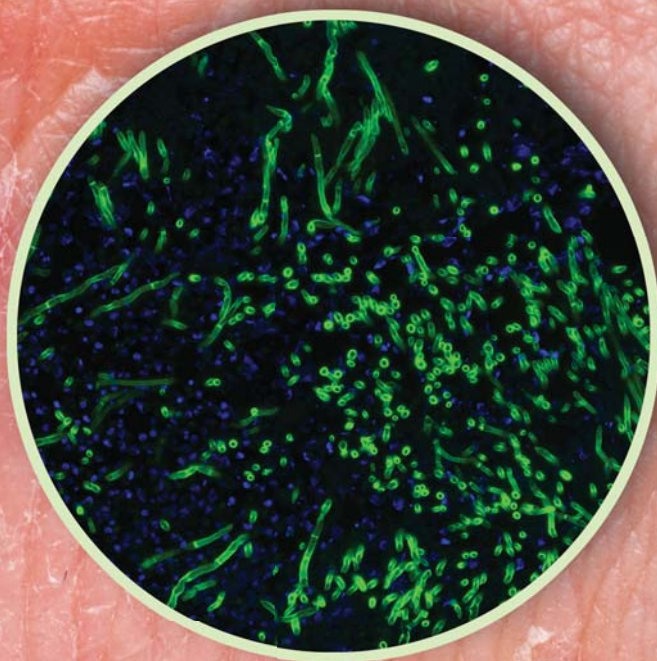

Fluorescence Detection of fungi in tissue

New Detection Method
of Pathogenic Fungi in
Tissuep2

Cronobacter spp.p3

Photo Competitionp6

Identifying Micro-
organisms Using Silver
Staining Techniquesp 7

Fast and Highly Specific Histological Detection of Pathogenic Fungi in Human Tissue

Monika Baeumle, PhD, Product Manager Biochemistry ...monika.baeumle@sial.com

Bernhard Schoenenberger, PhD, Supervisor R&D

Jakob Zbaeren, Thrombose Laboratory, Inselspital Bern

Lectins are ubiquitous proteins or glycoproteins that can be isolated from plant and animal sources and can bind to specific carbohydrate moieties.

Due to their high affinity to sugar residues, lectins have become important tools for sensitive detection of cellular carbohydrates, revealing subtle alteration in glycosylation between otherwise indistinguishable cells. This allows identification of cellular surface structures, e.g. cell surface, cytoplasm, and nuclear structures. Furthermore, lectin affinity binding allows for the detection of pathogenic degeneration of tissue as well as pathogenic infestations such as fungi.

Histochemical studies are of importance in the histological and pathological investigation of tissue in clinical research. Lectin histochemistry can be performed on living cells in suspension, on cell smears, tissue imprints, fixed tissue sections or fresh cryostat sections.

The recently developed Atto-dye labeled lectins have many applications, including carbohydrate, mitogenic and histochemical studies. Atto-dyes have very bright fluorescent signals and high photo stability, which enable a direct one step tissue-binding protocol. Time-consuming multistage amplification procedures are not required for Atto-dye lectin conjugates. Here, we demonstrate a highly specific identification of pathogenic fungi on human tissue via direct fluorescence detection using fluorescently labeled lectin (Figure 1).

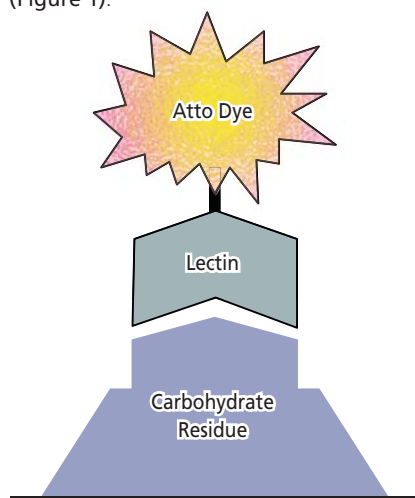


Figure 1: Direct one step binding of fluorescent labeled lectins.

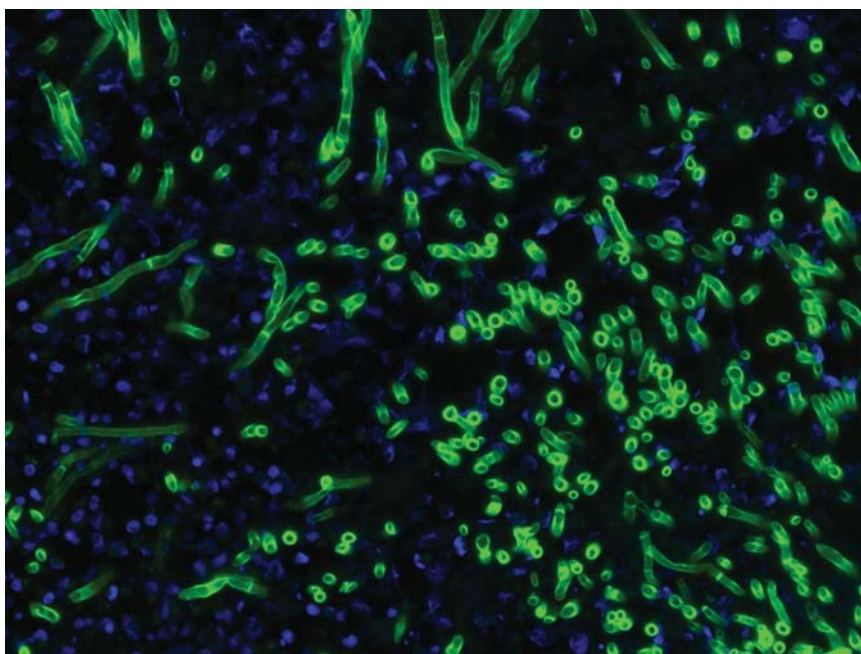


Figure 2: Fluorescence microscopy of human skin tissue section (paraffin fixation) with fungal infection. The target carbohydrate subunit chitotriose [(GlcNAc)₃] of the pathogenic fungi are specifically bound to lectin from *Phytolacca americana* Atto 488 conjugate (green). The nuclei are counterstained with DAPI (blue). Image by J. Zbaeren, Inselspital Bern, Switzerland.

Lectin histology was performed on both polymer and paraffin embedded human skin tissue. The lectin conjugate used was *Phytolacca americana*- Atto 488 (Cat. No. 39905). The conjugate was diluted 100 times in PBS buffer (pH 7.4) before incubating with each specimen for 30 min. After washing to remove any unbound lectin and counterstaining the nuclei with DAPI (Cat. No. 32670), the samples were examined using a microscope equipped for epifluorescence with a 450–490 nm excitation bandpass filter and a 520–560 nm barrier (emission) filter.

The images obtained show a very specific labeling of pathogenic fungi infecting human tissue (see Figure 2). The image demonstrates the fine filaments of the fungi containing typical *mycelium*, and individual fungi cells are clearly visible. A slightly higher fluorescence

is observed in the separating cross-walls between two cells (septa), which are due to a higher concentration of target carbohydrates. Very low background is observed.

Fungal cell walls contain chitin, a polymer of β -(1→4) linked *N*-acetyl-D-glucosamine, while animal and plant cells do not synthesize chitin. The lectin *Phytolacca americana* targets the fungal carbohydrate fragment chitotriose [(β -*N*-Acetyl-D-glucosamine)₃, (GlcNAc)₃] shown in green (λ_{ex} 485 nm). Due to the lack of the target carbohydrate chitotriose in the skin tissue, no specific interaction between the lectin *Phytolacca americana* and the tissue is observed. The bright and stable fluorescence properties of the Atto 488 dye provide a strong fluorescent signal without requiring additional amplification steps. Further experiments with staining different fungal in-

fectured tissues were carried out. Similar results confirm this approach to be a successful and reliable way to detect fungi. This application may encourage scientists to investigate further histological phenomena by using lectin interactions.

The Atto-dye lectin conjugates below are now available from Sigma Life Science. Additional lectins and lectin conjugates from Sigma Life Science may be found on at sigma-aldrich.com/enzymeexplorer.

Description	lex/lem (nm)	Carbohydrate Specificity	Cat. No.	Package Size
Concanavalin A - Atto 565-conjugate	563 / 592 in PBS	α -Mannose, α -Glucose	69535	1 mg
Lectin from Artocarpus integrifolia -Atto 594 conjugate	601/ 632 in PBS	O-Methyl- α -Galactose	76158	1 mg
Lectin from Ulex europaeus - Atto 488 conjugate	501 / 523 in PBS	α -L-Fucose	19337	0.5 mg
Lectin from Ulex europaeus - Atto 550 conjugate	554 / 576 in PBS	α -L-Fucose	94165	0.5 mg
Lectin from Ulex europaeus - Atto 594 conjugate	601/ 632 in PBS	α -L-Fucose	73873	0.5 mg
Lectin from Phaseolus vulgaris - Atto 488 conjugate (Leucoagglutinin)	501 / 523 in PBS	GlcNAc - Man	75319	1 mg
Lectin from Phaseolus vulgaris - Atto 550 conjugate (Leucoagglutinin)	554 / 576 in PBS	GlcNAc - Man	90852	1 mg
Lectin from Phaseolus vulgaris - Atto 647N conjugate (Leucoagglutinin)	644 / 669 in PBS	GlcNAc - Man	77363	1 mg
Lectin from Phytolacca Americana - Atto 488 conjugate	501 / 523 in PBS	(GlcNAc) ₃	39905	1 mg
Lectin from Phytolacca Americana - Atto 550 conjugate	554 / 576 in PBS	(GlcNAc) ₃	94816	1 mg
Lectin from Phytolacca Americana - Atto 647N conjugate	644 / 669 in PBS	(GlcNAc) ₃	03065	1 mg
Lectin from Triticum vulgaris - Atto 488 conjugate	501 / 523 in PBS	(GlcNAc) ₂ , α -N-acetylneuraminic acid	16441	1 mg
Lectin from Triticum vulgaris - Atto 532 conjugate	532 / 558 in PBS	(GlcNAc) ₂ , α -N-acetylneuraminic acid	68917	1 mg

Table 1: Fluorescence labeled lectins


Reference: Lectin Methods and Protocols; J. M. Rhodes; J. D. Milton; Humana Press, Totowa, New Jersey, 1997

Cronobacter spp. Classic and New Detection Methods

Jvo Siegrist, Product Manager Microbiology.... jvo.siegrist@sial.com

Enterobacter sakazakii, now reclassified as a novel genus called *Cronobacter*, is known to be related to neonatal infections.

Cronobacter is a rod-shaped, motile and facultatively-anaerobic bacteria of the family Enterobacteriaceae. Originally *Cronobacter* spp. were listed as yellow-pigmented *Enterobacter cloacae* (see **Table 1**). The bacteria was then later on called *Enterobacter sakazakii*. Taxonomic studies have determined that *E. sakazakii* comprises a high genetic heterogeneity and should be reclassified as a novel genus, "Cronobacter" [1].

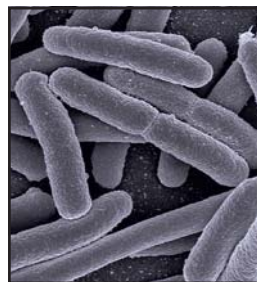

Kingdom: **Bacteria**
Phylum: **Proteobacteria**
Class: **Gamma Proteobacteria**
Order: **Enterobacteriales**
Family: **Enterobacteriaceae**
Genus: **Cronobacter**

Cronobacter spp. are ubiquitous and frequently found on vegetable, meat, fermented bread, dairy products and especially in baby food. Consumption of contaminated powdered infant formula milk (IFM) can result in sepsis, infant meningitis and necrotising enterocolitis. In most cases the victims are pre-term infants, low birth weight infants or immunocompromised infants in the first weeks. Mechanisms are not fully understood now but it is assumed that endotoxins are produced. The bacterium can adhere to and penetrate into various types of cells (e.g. endothelial cells). It also survives in macrophages.

As Enterobacteriaceae are susceptible to heat they do not survive most production processes and it is most likely that contamination happens after the process. The post process contamination can come from



Figure 1: Drinking baby

the addition of heat sensitive additives such as vitamins or other micronutrients or incorrect handling while reconstitution or storage. It has been shown, that stationary phase *Cronobacter* spp. are remarkably resistant to osmotic and dry stress compared with other species of the Enterobacteriaceae group. It is difficult to isolate *Cronobacter* spp. from samples as they are most likely stressed, unevenly distributed throughout the batch and also numbers will probably be fairly low, often lower than 1 CFU per g.

For the classical microbiological tests a pre-enrichment is used to recover the stressed cells followed by a selective enrichment step.

The FDA method recommends Enterobacteriaceae enrichment (EE or Mossel) broth which is then streaked onto VRBG agar and suspect colonies are subcultured onto TSA agar where the yellow pigmented colonies are confirmed by oxidase test and other biochemical tests. See also **Table 1** with all kind of biochemical reaction from *Cronobacter* spp. species (not only from FDA).

ISO/TS 22964:2008 methods recommends buffered peptone water (BPW) as pre-enrichment medium and modified lauryl sulphate broth with vancomycin (mLST) incubated at 44 °C for the secondary selective enrichment step. The next step is then a chromogenic agar for isolation and identification (see **Table 2**).

There are diverse chromogenic agars available which help to save work and time to result and are more reliable than traditional media. The detection principle is based on the alpha-glucosidase possessed by *Cronobacter* spp. (not by most other Enterobacteriaceae) which cleaves the 5-bromo-4-chloro-3-indolyl-a-D-glucopyranoside or similar substrates. The result is a plate with e.g. blue colonies in case of *Cronobacter* spp. (see **Figure 2** and **3**) but biochemical confirmation is still required.



Figure 2: HiCrome™ *Cronobacter* spp. Agar (Fluka 92324); *Cronobacter* spp. (blue), *E. aerogenes* (green) *K. pneumoniae* (yellow)



Figure 3: HiCrome™ *Cronobacter* spp. Agar, Modified (Fluka 14703), with *Cronobacter* spp. colonies

Today there are also studies that not all *Cronobacter* spp. giving yellow pigmented colonies on tryptic soy agar and it was showed that some type of strains did not grow at 44-45 °C. Also some selective media may be too selective to recover all the *Cronobacter* species. There is still some work to do to improve the official methods.



Did you know...

why this novel genus is called *Cronobacter*?

Cronos, was one of the Titans of the greek mythology who swallowed each of his children as soon as they were born. As *Cronobacter* species are harmful to neonates, the name was found to be adequate.

20% Test Discount
for *Cronobacter* spp.
media

(see table 2; Promo Code T62.
Valid until 30.09.2009)

Biochemical Test	Reaction of <i>Cronobacter</i>
Gram	-
oxidase	-
catalase	+
H ₂ S production	-
nitrate reduction	+
citrate utilization	+
esculin hydrolyzation	+
arginine hydolysation	+
Lysine	-
L-ornithine decarboxylation	+
Urease	-
Indole	-
ONPG	+
D-adonitol	-
L-arabinose	+
D-arabitol	-
D-cellobiose	+
Dulcitol	-/+
D-fructose	+
D-glucose	+
D-galactose	+
galacturonate	+
Inositol	+/-
Inulin	+
Lactose	+
Malonate	+/-
D-maltose	+
D-mannitol	+
D-mannose	+
D-melibiose	+
x-methyl-D-glucoside	+
D-raffinose	+
L-rhamnose	+
Salicin	+
Sorbitol	-
D-sucrose	+
D-trehalose	+
Xylose	+
Acetoin production (VP test)	+
methyl red test	-
tryptic soy agar at 25 °C	yellow pigmented

Table 1: Biochemical reactions from *Cronobacter* spp.

Brand	Cat. No.	Medium	Description
Fluka	92324	HiCrome™ <i>Cronobacter</i> spp. Agar	chromogenic media
Fluka	14703	HiCrome™ <i>Cronobacter</i> spp. Agar, Modified	chromogenic media according ISO
Fluka	69965	Mossel Broth (E.E. broth)	for pre-enrichment
Fluka	77187	Peptone Water, phosphate-buffered	for pre-enrichment
Fluka	22091	Tryptic Soy Agar	for isolation and differentiation
Fluka	79872	Tryptic Soy Agar	for isolation and differentiation

Table 2: Media for detection of *Cronobacter* spp.

A New Molecular Biology Method

Rapid detection and identification of *Cronobacter* species is required, since even low cell numbers have been reported to cause a disease. HybriScan® *D Cronobacter* spp. is a new rapid molecular test system for detection of bacteria of the genus *Cronobacter* in food, especially in dried infant formula milk and its production environment. It is based on the detection of rRNA by sandwich hybridisation and so **no PCR** is needed. It is a 96 well microplate format and the workflow is very similar to an ELISA test.

Figure 4 shows the specificity of HybriScan®- *Cronobacter* spp. Different cell amounts and related *Enterobacteriaceae* were tested within a validation study. No signals were observed using $2,3 \times 10^8$ *Enterobacter cloacae* cells or 7×10^8 *Citrobacter freundii* cells per assay, whereas clear specific signals were detectable using $2,6 \times 10^3$, $1,3 \times 10^4$, and $2,6 \times 10^4$ cells of *Cronobacter* species, respectively. These results demonstrate that the HybriScan® system is highly specific for *Cronobacter* spp.

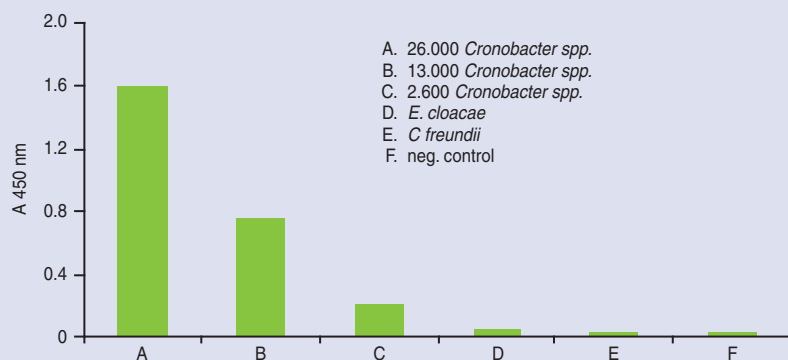


Figure 4: Specificity of HybriScan®- *Cronobacter* spp. Different cell numbers of *Cronobacter* spp. and related *Enterobacteriaceae* like *E. cloacae* and *Citrobacter freundii* were tested. Measurement data for HybriScan® analyses represent absorption at 450 nm.


A validation study of HybriScan®- *Cronobacter* spp. was performed using two different enrichment procedures: (1) single step enrichment for 24 – 26 hours at 37 °C in ESSB broth (*Enterobacter sakazakii* selective broth) and (2) two step enrichment starting with a pre-enrichment for 18 – 20 hours at 37 °C in buffered peptone water and followed by a selective enrichment for 24 – 26 hours at 45 °C in mLST selective broth. The results of the above mentioned validation study are presented in Figure 4.

Brand	Cat. No.	Name	Assays
Fluka	12838	HybriScan® <i>D Cronobacter</i> spp.	96

Table 2: Ordering Information

References:

1. Cawthorn, D.M.; Botha, S.; Witthuhn, R.S. Evaluation of different methods for the detection and identification of *Enterobacter sakazakii* isolated from South African infant formula milks and the processing environment. *International Journal of Food Microbiology*, 127:129–138 (2008)
2. Isolation and Enumeration of *Enterobacter sakazakii* from Dehydrated Powdered Infant Formula, U.S. Food and Drug Administration Center for Food Safety and Applied Nutrition (2002)
3. Color Atlas and Textbook of Diagnostic Microbiology, 5th edition, Lippincott Williams & Wilkins (1997)
4. Iversen et al., The taxonomy of *Enterobacter sakazakii*: proposal of a new genus *Cronobacter* gen. nov. and descriptions of *Cronobacter sakazakii* comb. nov., *Cronobacter sakazakii* subsp. *sakazakii*, comb. nov., *Cronobacter sakazakii* subsp. *malonicus* subsp. nov., *Cronobacter turicensis* sp. nov., *Cronobacter muytjensii* sp. nov., *Cronobacter dublinensis* sp. nov. and *Cronobacter* genomospecies 1, *BMC Evolutionary Biology*, 7:64 (2007)
5. Bergey's manual of determinative bacteriology, D.H. Bergey, J.G. Holt, 9th Edition, Lippincott Williams & Wilkins (1994)
6. M.B. Kleiman et al., Meningoencephalitis and Compartmentalization of the Cerebral Ventricles Caused by *Enterobacter sakazakii*, *Journal of Clinical Microbiology*, p. 352-354 (1981)
7. ISO/TS 22964:2006 Milk and milk products detection of *Enterobacter Sakazakii*
8. O. Guillaume-Gentil, V. Sonnard, M.C. Kandahai, J.D. Mauragg, H. Jootsen, A simple and Rapid Cultural Method for Detection of *Enterobacter Sakazakii* in environmental samples. *Journal of Food Protection*, 68 (1), 2005, pp. 64-69
9. K. Riedel, A. Lehner, Identification of proteins involved in osmotic stress response in *Enterobacter sakazakii* by proteomics. *Proteomics* 7, 1217-1231 (2007)
10. F.J. Pagotto, M. Nazarowec, S. Bidawid, J.M. Farber, *Enterobacter sakazakii*: Infectivity and enterotoxin production in vitro and in vivo; *J. of Food Protection*, Vol. 66, 3, p. 370-375 (2003)

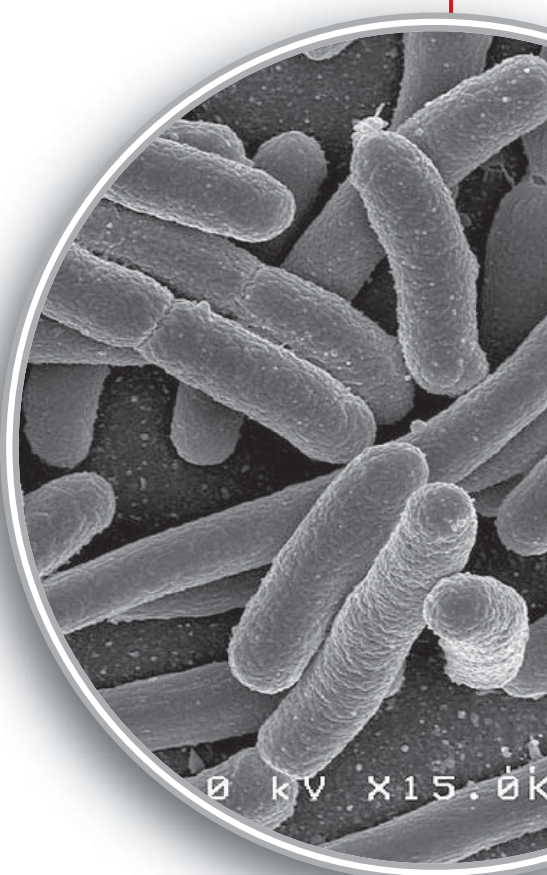


FLUKA-Microbiology Photography Competition!

General Information

This competition for photography is sponsored by Sigma-Aldrich with the aim of encouraging microbiologist to show something about their work and science.

The four best photographic entries will be presented in the Microbiology Focus and the *Best of Show* will get a place on a cover! All 4 winners will be awarded a trendy USB stick.

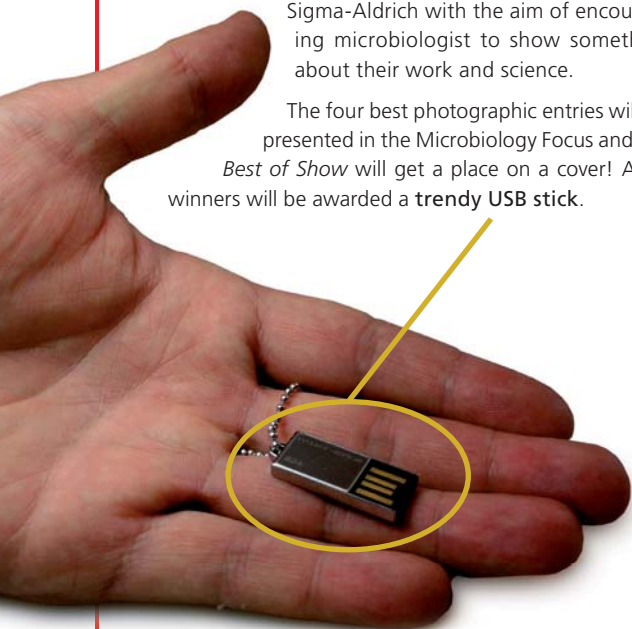

The competition will be judged by:

Prof. Mohammad Manafi
Medical University of Vienna,
Head of Department for Food Hygiene

Dr. Antje Breitenstein
Scanbec GmbH, CEO

Prof. Dr. Corinne Gantenbein
ZHSW, Head of Microbiology Department

Jvo Siegrist
Sigma-Aldrich, Product Manager Microbiology

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
Method of Entry

There is no entry fee. An individual may enter a maximum of 3 times and an entry form must be completed for each entry. To submit a photo, go to the following website:

sigma-aldrich.com/mibi_competition

Rules of the Competition and Conditions of Entry

1. The competition is open to all residents worldwide.
2. Entries should illustrate any microorganisms (living or dead) or a microbiologist in action at work
3. Picture size should be at least 400 dpi and 90 x 120 mm (max 3 MB). The file format must be in jpg, tiff or pdf.
4. The entries will be judged on:
 - clarity of presentation
 - composition
 - illumination and contrast
 - congruency of subject matter and title of photograph
 - scientific interest and relevance
 - originality
5. Winning entries will be retained by Sigma-Aldrich, who will have sole rights of publication, reproduction and display.
6. Closing date will be 30th Sept. 2009.
7. Entries after the closing date will not be considered. Entries received incomplete, illegible, mutilated, altered or not complying exactly with the instructions and theme may be disqualified.
8. Decisions of the judges in all matters affecting the competition will be final and legally binding.



sigma-aldrich.com/mibi_competition

Identifying Microorganisms Using Silver Staining Techniques

Mark Frei Technical Marketing Specialist.... mark.frei@sial.com

Silver staining techniques rely on basic chemical reactions for the microscopic examination and identification of microorganisms. These stains aid pathologists in the evaluation of disease states and help guide physicians in patient treatment.

Silver staining methods can be 10 – 100 times more sensitive than other staining techniques that rely on dyes that must penetrate the cell, however silver stains have a tendency to be capricious. Non-specific silver deposition and over-staining can produce a loss of detail. Clean glassware, proper technique and high quality reagents are necessary to obtain satisfactory results

Sigma-Aldrich supplies two modified silver staining kits that allow more reliable and consistent staining compared to the traditional method.

The Silver Stain Kit (Modified Steiner-Steiner) is intended for the demonstration of spirochetes and non-filamentous bacteria in sections of paraffin-embedded tissue. The method provides the option to utilise a microwave oven to accelerate and accentuate the silver stain in tissue sections. The heat produced in the microwave facilitates the impregnation of silver nitrate into the tissues, resulting in a much cleaner background than the traditional method. The organisms stain black and the background is yellow-brown (see **Figure 1**). Try the Modified Steiner-Steiner Kit (HT101-A) for consistent and reproducible staining results.

The Silver Stain Kit (modified GMS) is intended for use in histological visualisation of fungi, basement membrane and some opportunistic organisms. The organisms are stained black whilst other tissue elements are stained green (see **Figure 2**). Gomori Methenamine Silver (GMS) traditionally requires elaborate solution preparation and results can vary considerably due to the capricious nature of metal impregnation and photographic development. The Modified GMS Kit (HT100-A) incorporates a stable working silver methenamine salt along with buffer, toning reagent and developer, avoiding these issues. The use of a microwave oven allows for more rapid staining. Silver stains are of great importance in biomedical research applications and diagnostic pathology. Due to their low cost and ease of interpretation, silver stains will remain an important complement to emerging molecular biology technologies.

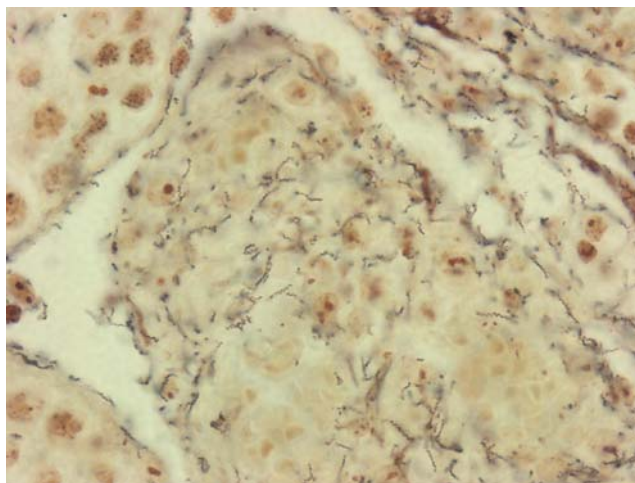


Figure 1: Steiner-Steiner silver stain of spirochetes


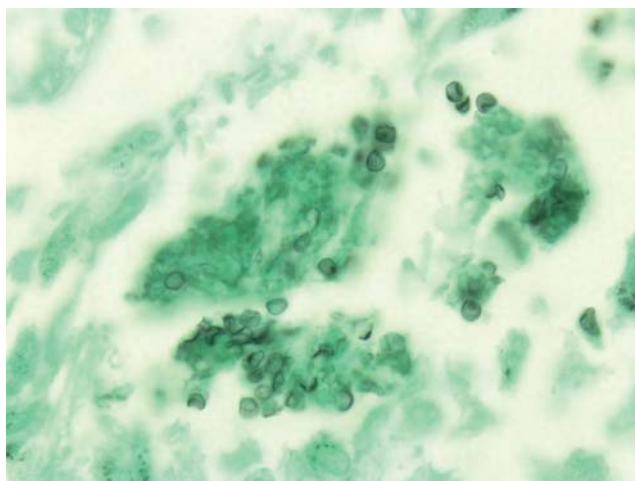


Figure 2: GMS silver stain of Pneumocystis

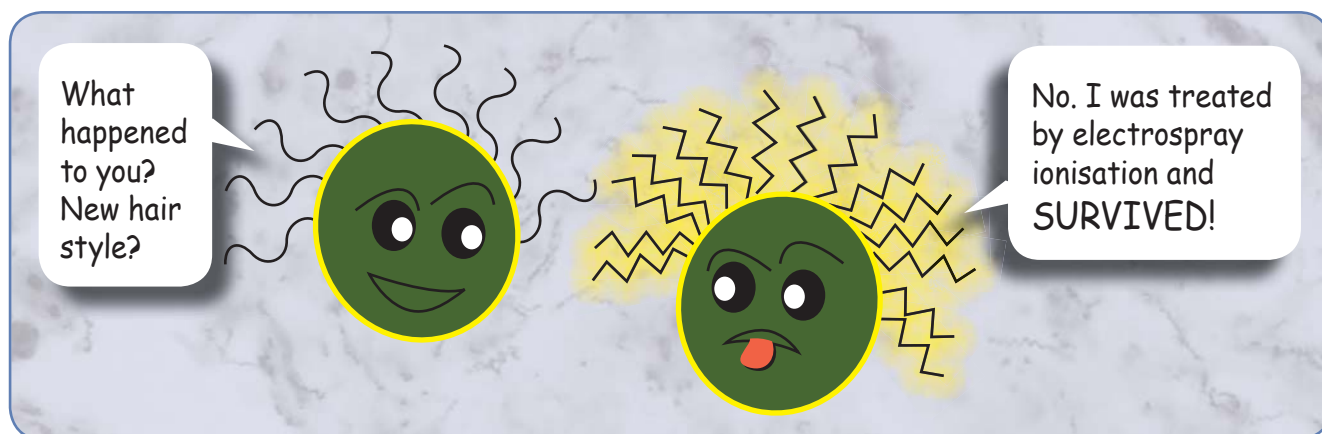

What bacteria think about new methods

Argentina

SIGMA-ALDRICH DE ARGENTINA S.A.
Free Tel: 0810 888 7446
Tel: (+54) 11 4556 1472
Fax: (+54) 11 4552 1698

Australia

SIGMA-ALDRICH PTY LTD.
Free Tel: 1800 800 097
Free Fax: 1800 800 096
Tel: (+61) 2 9841 0555
Fax: (+61) 2 9841 0500

Austria

SIGMA-ALDRICH HANDELS GmbH
Tel: (+43) 1 605 81 10
Fax: (+43) 1 605 81 20

Belgium

SIGMA-ALDRICH NV/S.A.
Free Tel: 0800 14747
Free Fax: 0800 14745
Tel: (+32) 3 899 13 01
Fax: (+32) 3 899 13 11

Brazil

SIGMA-ALDRICH BRASIL LTDA.
Free Tel: 0800 701 7425
Tel: (+55) 11 3732 3100
Fax: (+55) 11 5522 9895

Canada

SIGMA-ALDRICH CANADA LTD.
Free Tel: 1800 565 1400
Free Fax: 1800 265 3858
Tel: (+1) 905 829 9500
Fax: (+1) 905 829 9292

Chile

SIGMA-ALDRICH QUIMICA LIMITADA
Tel: (+56) 2 495 7395
Fax: (+56) 2 495 7396

China

SIGMA-ALDRICH (SHANGHAI)
TRADING CO. LTD.
Free Tel: 800 819 3336
Tel: (+86) 21 6141 5566
Fax: (+86) 21 6141 5567

Czech Republic

SIGMA-ALDRICH spol. s r. o.
Tel: (+420) 246 003 200
Fax: (+420) 246 003 291

Denmark

SIGMA-ALDRICH DENMARK A/S
Tel: (+45) 43 56 59 00
Fax: (+45) 43 56 59 05

Finland

SIGMA-ALDRICH FINLAND OY
Tel: (+358) 9 350 9250
Fax: (+358) 9 350 92555

France

SIGMA-ALDRICH CHIMIE S.à.r.l.
Free Tel: 0800 211 408
Free Fax: 0800 031 052
Tel: (+33) 474 82 28 00
Fax: (+33) 474 95 68 08

Germany

SIGMA-ALDRICH CHEMIE GmbH
Free Tel: 0800 51 55 000
Free Fax: 0800 64 90 000
Tel: (+49) 89 6513 0
Fax: (+49) 89 6513 1160

Greece

SIGMA-ALDRICH (O.M.) LTD.
Tel: (+30) 210 994 8010
Fax: (+30) 210 994 3831

Hungary

SIGMA-ALDRICH Kft
Ingyenes telefonszám: 06 80 355 355
Ingyenes fax szám: 06 80 344 344
Tel: (+36) 1 235 9055
Fax: (+36) 1 235 9050

India

SIGMA-ALDRICH CHEMICALS
PRIVATE LIMITED
Telephone
Bangalore: (+91) 80 6621 9400
New Delhi: (+91) 11 4358 8000
Mumbai: (+91) 22 2570 2364
Hyderabad: (+91) 40 4015 5488
Kolkata: (+91) 33 4013 8003
Fax
Bangalore: (+91) 80 6621 9650
New Delhi: (+91) 11 4358 8001
Mumbai: (+91) 22 2579 7589
Hyderabad: (+91) 40 4015 5466
Kolkata: (+91) 33 4013 8016

Ireland

SIGMA-ALDRICH IRELAND LTD.
Free Tel: 1800 200 888
Free Fax: 1800 600 222
Tel: (+353) 402 20370
Fax: (+353) 402 20375

Israel

SIGMA-ALDRICH ISRAEL LTD.
Free Tel: 1 800 70 2222
Tel: (+972) 8 948 4100
Fax: (+972) 8 948 4200

Italy

SIGMA-ALDRICH S.r.l.
Numero Verde: 800 827018
Tel: (+39) 02 3341 7310
Fax: (+39) 02 3801 0737

Japan

SIGMA-ALDRICH JAPAN K.K.
Tel: (+81) 3 5796 7300
Fax: (+81) 3 5796 7315

Korea

SIGMA-ALDRICH KOREA
Free Tel: (+82) 80 023 7111
Free Fax: (+82) 80 023 8111
Tel: (+82) 31 329 9000
Fax: (+82) 31 329 9090

Malaysia

SIGMA-ALDRICH (M) SDN. BHD
Tel: (+60) 3 5635 3321
Fax: (+60) 3 5635 4116

Mexico

SIGMA-ALDRICH QUÍMICA, S.A. de C.V.
Free Tel: 01 800 007 5300
Free Fax: 01 800 712 9920
Tel: (+52) 722 276 1600
Fax: (+52) 722 276 1601

The Netherlands

SIGMA-ALDRICH CHEMIE BV
Free Tel: 0800 022 9088
Free Fax: 0800 022 9089
Tel: (+31) 78 620 5411
Fax: (+31) 78 620 5421

New Zealand

SIGMA-ALDRICH NEW ZEALAND LTD.
Free Tel: 0800 936 666
Free Fax: 0800 937 777
Tel: (+61) 2 9841 0555
Fax: (+61) 2 9841 0500

Norway

SIGMA-ALDRICH NORWAY AS
Tel: (+47) 23 17 60 00
Fax: (+47) 23 17 60 10

Poland

SIGMA-ALDRICH Sp. z o.o.
Tel: (+48) 61 829 01 00
Fax: (+48) 61 829 01 20

Portugal

SIGMA-ALDRICH QUÍMICA, S.A.
Free Tel: 800 202 180
Free Fax: 800 202 178
Tel: (+351) 21 924 2555
Fax: (+351) 21 924 2610

Russia

SIGMA-ALDRICH RUS, LLC
Tel: (+7) 495 621 5579
Fax: (+7) 495 621 5923

Singapore

SIGMA-ALDRICH PTE. LTD.
Tel: (+65) 6779 1200
Fax: (+65) 6779 1822

Slovakia

SIGMA-ALDRICH spol. s r. o.
Tel: (+421) 255 571 562
Fax: (+421) 255 571 564

South Africa

SIGMA-ALDRICH (PTY) LTD.
Free Tel: 0800 1100 75
Free Fax: 0800 1100 79
Tel: (+27) 11 979 1188
Fax: (+27) 11 979 1119

Spain

SIGMA-ALDRICH QUÍMICA, S.A.
Free Tel: 900 101 376
Free Fax: 900 102 028
Tel: (+34) 91 661 99 77
Fax: (+34) 91 661 96 42

Sweden

SIGMA-ALDRICH SWEDEN AB
Tel: (+46) 8 742 4200
Fax: (+46) 8 742 4243

Switzerland

SIGMA-ALDRICH CHEMIE GmbH
Free Tel: 0800 80 00 80
Free Fax: 0800 80 00 81
Tel: (+41) 81 755 2828
Fax: (+41) 81 755 2815

United Kingdom

SIGMA-ALDRICH COMPANY LTD.
Free Tel: 0800 717 181
Free Fax: 0800 378 785
Tel: (+44) 1747 833 000
Fax: (+44) 1747 833 313

United States

SIGMA-ALDRICH
Toll-Free: 800 325 3010
Toll-Free Fax: 800 325 5052
Tel: (+1) 314 771 5765
Fax: (+1) 314 771 5757

Vietnam

SIGMA-ALDRICH PTE LTD. VN R.O.
Tel: (+84) 3516 2810
Fax: (+84) 6258 4238

Internet

sigma-aldrich.com



Mixed Sources
Product group from well-managed
forests, controlled sources and
recycled content fiber
www.fsc.org Cert no. SGS-COC-XXXXXX
© 1996 Forest Stewardship Council



World Headquarters

3050 Spruce St., St. Louis, MO 63103
(314) 771-5765
sigma-aldrich.com

Order/Customer Service (800) 325-3010 • Fax (800) 325-5052

Technical Service (800) 325-5832 • sigma-aldrich.com/techservice

Development/Custom Manufacturing Inquiries **SAFC®** (800) 244-1173

©2009 Sigma-Aldrich Co. All rights reserved. SIGMA, SAFC, **SAFC®**, SIGMA-ALDRICH, ALDRICH, FLUKA, SUPELCO, are trademarks belonging to Sigma-Aldrich Co. and its affiliate Sigma-Aldrich Biotechnology, L.P. Sigma brand products are sold through Sigma-Aldrich, Inc. Sigma-Aldrich, Inc. warrants that its products conform to the information contained in this and other Sigma-Aldrich publications. Purchaser must determine the suitability of the product(s) for their particular use. Additional terms and conditions may apply. Please see reverse side of the invoice or packing slip.

*Accelerating Customers'
Success through Innovation and
Leadership in Life Science,
High Technology and Service*